

## SUPPORT

0.65X


1X


0.5X

THE C-MOUNT - A NEEDFUL THING

**BASIC MICROSCOPY CONCEPTS** 



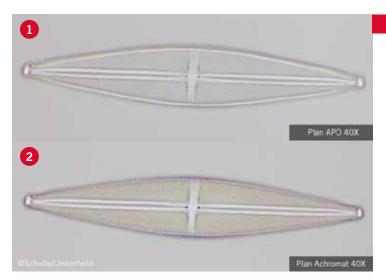

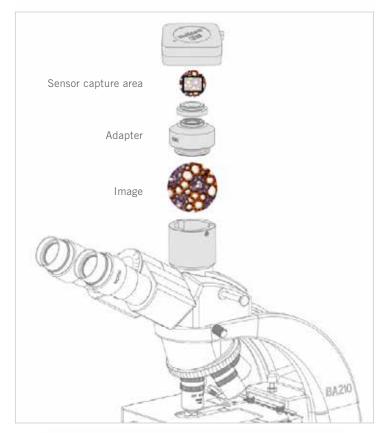



## THE C-MOUNT, A NEEDFUL THING

For beginners, the combination of a light microscope with a digital camera includes some surprises. There are obvious discrepancies between the visual impression in the eyepieces and on the screen respectively. These discrepancies are about color impression and image section.

The color impression in the eyepieces may be taken as a reference and depends on the color correction of the objective. The **Plan Apochromatic lens** design gives best preconditions for **maximal color fidelity and resolution**. The imperfection of a secondary spectrum, caused by the usage of white light, is compensated for 3 colors, while spherical aberrations are corrected for 2 wavelengths. The result **(1)** is an impressive image quality without any colored fringes. Especially in colorless black/white samples the absence of secondary spectrum becomes obvious. However, for price reasons, it is the **Plan Achromatic** quality level **(2)** of a modern Infinity System which sets the point of reference for most users.


Many beginners expect a 100% color transfer to the screen via digital camera. But the sensory hardware of the digital camera, a CMOS or CCD chip, acts as a bottleneck. CCD targets are superior to CMOS sensors in terms of color reproduction, but they still lack the dynamic range of the human eye. If it comes to the price point, a CMOS sensor will find his friends. A clever signal processing will lead to a reasonable image quality at an affordable price point.

There is one more problem to be solved, as the circular intermediate image within the microscope has to be captured by a rectangular sensor. Microscope manufacturers tend to enlarge the visible area (diameter of the intermediate image = Field Of View; FOV). Chip manufacturers go the opposite direction as the consumer market for pocket cameras and mobile phones requires small-sized sensors.

The combination of a typical microscope camera with the photo port of a microscope requires an intermediate piece of hardware, combining optical and mechanical tasks : the C-mount adapter (3).





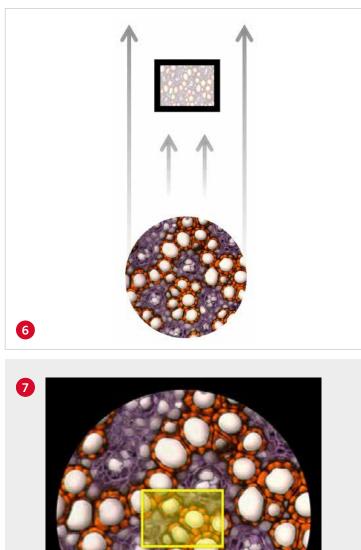

Adapter 1X - Captured area = 10% of FOV 20

Its upper part carries a standardized thread of 1 inch outer diameter and 1/32 inch pitch, ready for the mounting of the camera (4).

Its lower part refers to to the microscope manufacturer and the respective microscope model **(5)**. Despite some homogenisations, even within one brand there might be different interfaces for different microscopes.


A C-mount 1X without lens is a pure mechanical device for a direct positioning of the camera sensor "on top" of the intermediate image **(6)**.

This combination gives a disappointing image section of 10% Field of View (7).


A camera adapter with a magnification factor  $\leq 1$  will help. The adapter's optics "compresses" the intermediate image and thus increases the captured area (8).






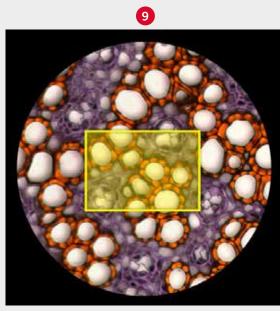

Adapter 0.65X - Captured area = 23% of FOV 20

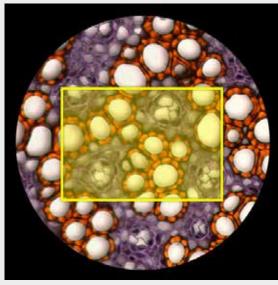


Adapter 0.5X - Captured area = 38% of FOV 20

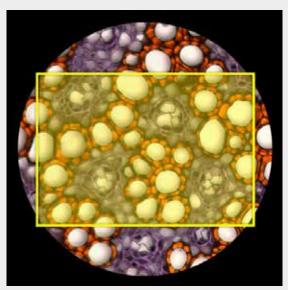


Adapter 0.35X - Captured area = 76% of FOV 20

Common "magnifications" are 0.65X, 0.5X and 0.35X. Compare the difference between the captured areas depending on the adapter chosen (9).

A C-mount with reducing lens simulates a larger sensor size, as shown in this comparison:

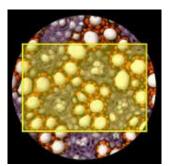

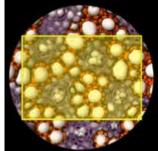

Sensor 1/2" Adapter 0.35X 76% FOV

Sensor 2/3" Adapter 0.5X 72% FOV

The disadvantage of a reducing C-mount should be taken into consideration. Any compression of the intermediate image reduces resolution as image structures are pushed together: With Great Compression comes Great Reduction (Uncle Ben to Peter Parker). Selecting the camera adapter, the user takes his decision for image section or resolution. That kind of dilemma is well known in microscopy: remember the setting of the aperture diaphragm which always is a compromise between contrast and resolution.



Canada | China | Germany | Spain | USA





If you want to know more about our products, visit our **Support Zone** at www.moticeurope.com/support

\*CCIS® is a trademark of Motic Incorporation Ltd. Motic Incorporation Limited Copyright © 2002-2016. All Rights Reserved.

Design Change: The manufacturer reserves the right to make changes in instrument design in accordance with scientific and mechanical progress, without notice and without obligation. Designed in Barcelona (Spain) August 2016

