





HOW TO MOUNT THE MOTICAM 1080 INT ONTO A BA SERIES MICROSCOPE



ideo version available in our Youtube channel. ww.voutube.com/MoticEurope








STEPS TO FOLLOW - Camera assembly

(1) (2)

Mount the screen to the camera and tighten the screw.

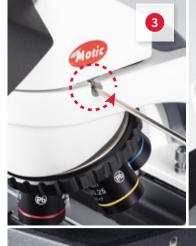



STEPS TO FOLLOW - Mount onto a microscope

(3) (4)

Loosen the screw with the allen key and remove the microscope head.

(5)

Attach the camera to the microscope body and tighten the screw.



(6)

Remove the black cover of the camera by loosening the screw.





(7)

Attach the microscope head on the camera by tightening the screw.

(8)

Turn the microscope and connect the HDMI cable to the HDMI ports of the camera and screen.

(9)

Connect the power cable coming from the camera into the DC-12V port of the screen.

(10)


10

Connect the power supply cable into the INPUT DC 12V port of the camera.



Canada | China | Germany | Spain | USA



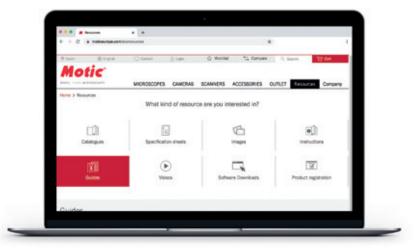

If you want to know more about our products, visit our **Resources Zone** at www.moticeurope.com/resources

*CCIS® is a trademark of Motic Incorporation Ltd. Motic Incorporation Limited Copyright © 2002-2021. All Rights Reserved.

Design Change: The manufacturer reserves the right to make changes in instrument design in accordance with scientific and mechanical progress, without notice and without obligation. Created and designed in Barcelona (Spain) by Motic Europe. February 2021.

