



BASIC MICROSCOPY CONCEPTS



SUPPORT

PARFOCALITY ADJUSTMENT FOR UPRIGHT MICROSCOPES





PARFOCALITY ADJUSTMENT

FOR UPRIGHT MICROSCOPES

Parfocality is a property of the optical components that allows the microscope to stay in focus when changing between lenses of different magnification.

This means that when the objectives of a microscope are changed from higher to lower magnification or vice versa, the sample you are viewing stays in focus.

This is important especially when taking photographs or making videos using a microscope and a camera. If the parfocality is not adjusted, the camera will not be focusing correctly, even if the image seen through the eyepieces is in focus.

To ensure the correct parfocality adjustment of your microscope, just follow these easy steps:

(1) Adjust the interpupillary distance so that both the right and left field of view become one.

(2) Set the diopter adjustment on both eyepieces (or on eyepieces tubes) on the "0" position.

(3) Select the lowest magnification objective and focus the image of the specimen with the help of the coarse/fine focusing knobs, using only one eye.* *Use the eye with which you're most comfortable with.

(4) Once you obtain the best possible focus, close the eye you just used and use your other eye for the next steps.

(5) Correct the focus for the second eye by using only the diopter adjustment.

Important: In this case do not use the coarse/fine focusing knobs.





(6) Now select the highest magnification and repeat the procedure to match the sharpness for higher magnification.

(7) Keep this final diopter position for all magnifications/ lenses. The diopter position of each user is different, for this reason you might want to take note of your graduation (from the scale), for easier resetting.

(8) The focus of the image should now be constant as you move to higher or lower magnification (parfocal).



Canada | China | Germany | Spain | USA





If you want to know more about our products, visit our **Support Zone** at www.moticeurope.com/support

*CCIS® is a trademark of Motic Incorporation Ltd. Motic Incorporation Limited Copyright © 2002-2016. All Rights Reserved.

Design Change: The manufacturer reserves the right to make changes in instrument design in accordance with scientific and mechanical progress, without notice and without obligation. Designed in Barcelona (Spain) November 2016

